Data Acquisition Toolbox™ SDK User's Guide

MATLAB®



R2017a

How to Contact MathWorks



The MathWorks, Inc. 3 Apple Hill Drive Natick, MA 01760-2098

Data Acquisition Toolbox[™] SDK User's Guide

© COPYRIGHT 2017 by The MathWorks, Inc.

The software described in this document is furnished under a license agreement. The software may be used or copied only under the terms of the license agreement. No part of this manual may be photocopied or reproduced in any form without prior written consent from The MathWorks, Inc.

FEDERAL ACQUISITION: This provision applies to all acquisitions of the Program and Documentation by, for, or through the federal government of the United States. By accepting delivery of the Program or Documentation, the government hereby agrees that this software or documentation qualifies as commercial computer software or commercial computer software documentation as such terms are used or defined in FAR 12.212, DFARS Part 227.72, and DFARS 252.227-7014. Accordingly, the terms and conditions of this Agreement and only those rights specified in this Agreement, shall pertain to and govern the use, modification, reproduction, release, performance, display, and disclosure of the Program and Documentation by the federal government (or other entity acquiring for or through the federal government) and shall supersede any conflicting contractual terms or conditions. If this License fails to meet the government's needs or is inconsistent in any respect with federal procurement law, the government agrees to return the Program and Documentation, unused, to The MathWorks, Inc.

Trademarks

MATLAB and Simulink are registered trademarks of The MathWorks, Inc. See www.mathworks.com/trademarks for a list of additional trademarks. Other product or brand names may be trademarks or registered trademarks of their respective holders.

Patents

MathWorks products are protected by one or more U.S. patents. Please see www.mathworks.com/patents for more information.

Revision History

March 2017 Online only

New for Version 3.11 (Release 2017a)



Contents

SDK Overview

Toolbox and Adaptor Hardware Driver Adaptor Hardware Driver Adaptor Hardware Driver Adaptor	
SDK Contents	1-3
Adaptor Creation Summary Tips	

Explore the SDK Demo Adaptor

2

1

Demo Adaptor Description	2-2
Source Files	2-2
Class Definitions in MATLAB	2-3
Executables	2-3
Enable the Demo Adaptor	2-5
Session Workflows with the Demo Adaptor	2-6
Device Discovery and Configuration	2-6
Single Scan Input and Output	2-8
Streaming Input and Output	2-10
Test the Demo Adaptor	2-12
Run Individual Tests	2-12
Run a Test Suite	2-13

Create Your Adaptor from the Demo Adaptor Edit and Build Your Adaptor	3-2 3-2
Use Your Adaptor in a Session	3-2 3-4
Modify Demo Tests for Your Adaptor	3-7
Further Suggestions	3-10
Errors and Exceptions	3-11
Nonstreaming	3 - 11
Streaming	3-11
Channel Groups	3-13
Channel Group Description	3 - 13
Channel Group Restrictions	3-14
Custom Functions	3-15
Vendor Adaptor Templates	3-17
Typical Workflow to Create Adaptor	3-17
Deliver Your Adaptor	3-19

API Reference

4

3

Adaptor API Reference
Lifetime
Enumeration
Hardware Management
Vendor and Device Discovery
Subsystem Discovery
Configuration
Reservation
SingleScans

Streaming API Reference	4-21
Initialization and Configuration	4-21
Start and Stop	4-22
Data Availability	4-23
Transfer Data	4-2 4

State and Sequence Diagrams

5

6

State Machine Diagram		
Streaming Sequence Diagrams	5-4	
Foreground Streaming Sequences	5-5	
Sequence for Finite Foreground Input	5-5	
Sequence for Finite Foreground Output	5-7	
Sequence for Finite Foreground Duplex Channel	5-9	
Background Streaming Sequences	5-12	
Sequence for Finite Background Input	5 - 12	
Sequence for Continuous Background Input with Stop	5 - 14	
Sequence for Finite Background Input with Wait	5-16	
Sequence for Finite Background Input with Stop Race	5-18	
Sequence for Errors and Exceptions	5-21	

Functions — Alphabetical List


SDK Overview

- "Toolbox and Adaptor" on page 1-2
- "SDK Contents" on page 1-3
- "Adaptor Creation Summary" on page 1-4

Toolbox and Adaptor

Hardware Driver Adaptor

The hardware driver adaptor is the interface between the data acquisition engine and the hardware driver. The adaptor's main purpose is to pass information between $MATLAB^{\circledast}$ and your hardware device via its driver.

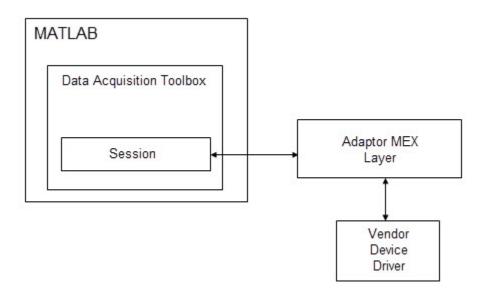

Hardware drivers are provided by the device vendor. For example, to acquire data using a National Instruments[®] board, the appropriate version of the driver must be installed on your system. Hardware drivers are not installed as part of the toolbox, but a suitable driver is usually installed on PCs that are equipped with a sound card. For any other devices, the drivers must be installed.

Related Examples

• "Adaptor Creation Summary" on page 1-4

SDK Contents

An installation of Data Acquisition ToolboxTM includes the following folders to support its SDK. Your MATLAB installation location is referred to as *matlabroot*.

Folder	Description
<pre>matlabroot\toolbox\daq\daqsdk\bin \win64</pre>	Built executable files for demo and vendor adaptors.
<pre>matlabroot\toolbox\daq\daqsdk \tests\+daq\+sdk\+tests</pre>	Test files for adaptors. The installed set of test files is for the demo adaptor.
<i>matlabroot</i> \toolbox\daq\daqsdk\src \daqadaptor	Adaptor C++ source files, one folder for each adaptor, and one other folder named Shared for elements common to all adaptors. DemoAdaptor contains all the source files for the demo adaptor. VendorAdaptor contains a set of templates.

Related Examples

- "Adaptor Creation Summary" on page 1-4
- "Create Your Adaptor from the Demo Adaptor" on page 3-2
- "Modify Demo Tests for Your Adaptor" on page 3-7
- "Vendor Adaptor Templates" on page 3-17

Adaptor Creation Summary

This topic provides a summary of adaptor creation with the SDK. For details and examples of these steps, see "Create Your Adaptor from the Demo Adaptor" on page 3-2.

- 1 Copy the demo adaptor or vendor adaptor source files into your working folder.
- 2 Change the names of the source files to reflect your own adaptor name.
- **3** Update the content of the source files so that the new names are used for references to other files, the adaptor, devices, and vendor.
- **4** Update the source file functions to use your driver code. For more information, see "Adaptor API Reference" on page 4-2.
- **5** Build the adaptor with the buildAdaptor function. Add the folder containing the built MEX-file to your MATLAB path.
- **6** Copy the demo adaptor tests and modify them for your adaptor. Add the test package folder to your MATLAB path.
- 7 Run the tests on your adaptor.
- 8 Deliver the finished adaptor MEX-file with your device driver and supporting files.

Tips

- Update, build, and test your adaptor iteratively one step at a time. Develop and test in small increments, proceeding upon the success of each step.
- When modifying the source files, do not remove any of the functions. Even if you do not use all the functions, they must be present when using buildAdaptor.

Related Examples

- "Create Your Adaptor from the Demo Adaptor" on page 3-2
- "Modify Demo Tests for Your Adaptor" on page 3-7

More About

- "Vendor Adaptor Templates" on page 3-17
- "Deliver Your Adaptor" on page 3-19

Explore the SDK Demo Adaptor

- "Demo Adaptor Description" on page 2-2
- "Enable the Demo Adaptor" on page 2-5
- "Session Workflows with the Demo Adaptor" on page 2-6
- "Test the Demo Adaptor" on page 2-12

Demo Adaptor Description

The demo adaptor installed with Data Acquisition Toolbox consists of the files described in the following tables.

In this section...

"Source Files" on page 2-2

"Class Definitions in MATLAB" on page 2-3

"Executables" on page 2-3

Source Files

The demo adaptor source files are in *matlabroot*\toolbox\daq\daqsdk\src \daqadaptor\DemoAdaptor.

File	Description
demoadaptor.hpp,demoadaptor.cpp	Wraps device driver code in methods that allow you to configure, discover, and enumerate the hardware in MATLAB.
daqstream_analog.hpp, daqstream_analog.cpp	Implements a DAQStream object for an analog subsystem, that allows you to stream data to and from the hardware.
custom_demo.cpp	Dispatches calls from MATLAB to custom functions in the demo adaptor. At a minimum this must contain a customizeMap function.

In addition to these files, the demo adaptor also uses some of the source files in *matlabroot*\toolbox\daq\daqsdk\src\daqadaptor\Shared:

Files	Purpose
adaptorfactory.cpp, adaptorfactory.hpp	Create adaptor for dispatch and streaming.
daqadaptor.cpp, daqadaptor.hpp	Implement adaptor class.
daqapi.h	C interface.

Files	Purpose
daqdatatypes.hpp	C++ equivalents of session data types.
daqinterfaces.hpp	IAdaptor/IDriver, called before streaming.
daqstream.cpp, daqstream.hpp	Transfer streaming data between MATLAB and device driver
dispatcher.cpp,dispatcher.hpp, dispatcher_common.hpp	MATLAB calls to convert data and call adaptor functions.
fakevendordriver.hpp	Fake or virtual driver for testing and demonstrations.
globals.h	Global settings.
mxconvert.hpp	Utility functions for data type conversions.

Class Definitions in MATLAB

The demo adaptor class definitions are in *matlabroot*\toolbox\daq\daqsdk\+daq\+demoadaptor.

File	Description
Session.m	Defines daq.demoadaptor.Session class.
	Defines vendor driver class for daq.getVendor.

In addition to the files in this table, the demo adaptor also uses some of the class definition files in *matlabroot*\toolbox\daq\daqsdk\+daq\+sdk.

Executables

The following demo adaptor executables are in *matlabroot*\toolbox\daq\daqsdk \bin\win64.

File	Description
DemoAdaptor.mexw64	Built demo adaptor.
	Accommodates streaming channel communication.

File	Description
daqmlconverter.dll	Handles data type conversion.

See Also

Functions enableDemoAdaptorDiscovery

Related Examples

- "Enable the Demo Adaptor" on page 2-5
- "Session Workflows with the Demo Adaptor" on page 2-6
- "Test the Demo Adaptor" on page 2-12

Enable the Demo Adaptor

By default, the demo adaptor is disabled when first installed. Use the following MATLAB command to enable it.

daq.sdk.utility.enableDemoAdaptorDiscovery

The adaptor is now ready for use. Confirm this with the command:

daq.getVendors

The output includes an entry for the demo adaptor with the vendor ID of mw:

index ID Operational Comment
1 mw true MathWorks

See Also

Functions enableDemoAdaptorDiscovery

Related Examples

- "Demo Adaptor Description" on page 2-2
- "Session Workflows with the Demo Adaptor" on page 2-6
- "Test the Demo Adaptor" on page 2-12

Session Workflows with the Demo Adaptor

In this section...

"Device Discovery and Configuration" on page 2-6 "Single Scan Input and Output" on page 2-8 "Streaming Input and Output" on page 2-10

Device Discovery and Configuration

When you create a data acquisition session, it applies to a specific vendor, and allows you to add applicable devices and channels. Discovery and configuration is part of setting up your session. This example shows a typical session setup with the demo adaptor.

Note To enable the demo adaptor in your installation, see the instructions in "Enable the Demo Adaptor" on page 2-5.

```
v = daq.getVendors
v =
Number of vendors: 2
index ID Operational Comment
1 ni false Click here for more info
2 mw true MathWorks
```

Properties, Methods, Events

Additional data acquisition vendors may be available as downloadable support packages. Open the Support Package Installer to install additional vendors.

```
d = daq.getDevices
```

d =

Data acquisition devices:

index Vendor Device ID Description

1	mw	MWDev0	MathWorks	MW314159
2	mw	MWDev1	MathWorks	MW314159
3	mw	MWDev2	MathWorks	MW628318

With a listing of available vendors and devices, you can create a session and add channels to it.

```
s = daq.createSession('mw')
```

s =

```
Data acquisition session using MathWorks hardware:
Will run for 1 second (1000 scans) at 1000 scans/second.
No channels have been added.
```

Properties, Methods, Events

To see details about one of the devices, use its index to access the array of devices.

d(1)

```
mw: MathWorks MW314159 (Device ID: 'MWDev0')
Analog input subsystem supports:
    3 ranges supported
    Rates from 0.1 to 1000000.0 scans/sec
    2 channels ('ai0','ai1')
    'Voltage','Current' measurement types
Analog output subsystem supports:
    3 ranges supported
    Rates from 0.1 to 1000000.0 scans/sec
    2 channels ('ao0','ao1')
    'Voltage','Current' measurement types
```

The first and second devices are the same model, so this example uses one for input ('ai0' and the other for output ('a00').

```
ch1 = addAnalogInputChannel(s,'MWDev0','ai0','voltage')
ch1 =
Data acquisition analog input voltage channel 'ai0' on device 'MWDev0':
```

```
Coupling: DC
 TerminalConfig: Differential
          Range: -10 to +10 Volts
           Name: ''
             ID: 'ai0'
         Device: [1x1 dag.sdk.DeviceInfo]
MeasurementType: 'Voltage'
ch2 = addAnalogOutputChannel(s, 'MWDev1', 'ao0', 'voltage')
ch2 =
Data acquisition analog output voltage channel 'ao0' on device 'MWDev1':
 TerminalConfig: SingleEnded
          Range: -10 to +10 Volts
           Name: ''
             ID: 'ao0'
         Device: [1x1 dag.sdk.DeviceInfo]
MeasurementType: 'Voltage'
```


View the session to see the configuration.

```
s
s =
Data acquisition session using MathWorks hardware:
  No data queued. Will run at 1000 scans/second.
  Number of channels: 2
     index Type Device Channel MeasurementType
                                                        Range
                                                                   Name
      1
           ai
              MWDev0 ai0
                               Voltage (Diff)
                                                  -10 to +10 Volts
     2
           ao MWDev1 ao0
                               Voltage (SingleEnd) -10 to +10 Volts
```

The session is now ready to send and receive single scans or streams of data.

Single Scan Input and Output

A single scan is when you send an output or read input from the channels at one moment in time. The data transfer is handled by the adaptor MEX layer. For the demo adaptor this is contained in the MEX-file matlabroot\toolbox\daq\daqsdk\bin \win64\DemoAdaptor.mexw64, which provides the functionality shown in the following diagram.

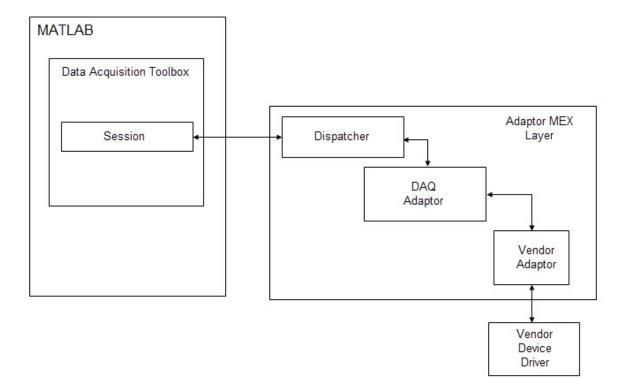

To generate a single scan analog output of 1.25 V, enter the following code.

```
outputSingleScan(s,1.25)
```

DemoDriver output: 1.25

To read a single scan of analog input:

```
data = inputSingleScan(s)
```

With only one input channel, this returns only a single value. If you add more analog input channels to the session, inputSingleScan returns a vector, with an element for each channel.

```
ch3 = addAnalogInputChannel(s,'MWDev2','ai0','voltage');
data = inputSingleScan(s)
data =
```


0 200

Remove channels from the session you no longer need.

removeChannel(s,[1 2 3]);

Streaming Input and Output

Streaming involves a sequence of input or output data on each channel, typically a waveform, comprised of many scans. Streaming can be accomplished in the foreground (blocking MATLAB until the stream is complete), or in the background (running asynchronously while MATLAB continues). Streaming might involve more scans or samples than the device memory can hold. For these reasons, the toolbox uses streaming channels to accommodate data flow. This allows data to be sent and received without causing a memory overflow, and without interrupting MATLAB.

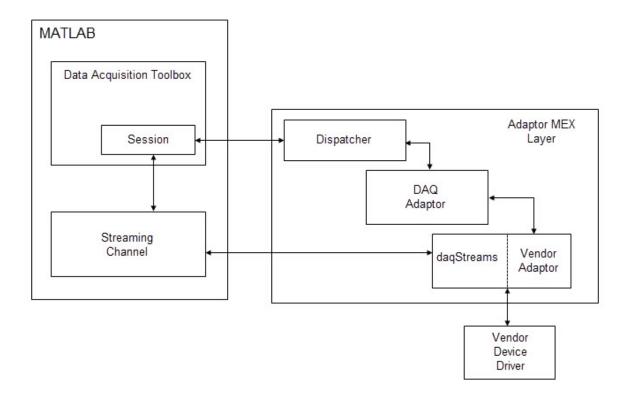

Use MWDev2 to generate a 1 KHz sinewave in the background for ten seconds. The default sample rate is 1000 per second.

```
ch4 = addAnalogOutputChannel(s, 'MWDev2', 'ao1', 'voltage')
ch4 =
Data acquisition analog output voltage channel 'ao1' on device 'MWDev2':
 TerminalConfig: SingleEnded
          Range: -10 to +10 Volts
           Name: ''
             ID: 'ao1'
         Device: [1x1 daq.sdk.DeviceInfo]
MeasurementType: 'Voltage'
Y = sin(linspace(0,2*pi*10,10000))' % 10 cycles in 10000 samples;
queueOutputData(s,Y);
startBackground(s)
s.IsRunning
  logical
   1
pause (10)
s.IsRunning
  logical
   0
```

See Also

Functions enableDemoAdaptorDiscovery

Related Examples

- "Demo Adaptor Description" on page 2-2
- "Enable the Demo Adaptor" on page 2-5
- "Test the Demo Adaptor" on page 2-12

Test the Demo Adaptor

In this section ...

"Run Individual Tests" on page 2-12

"Run a Test Suite" on page 2-13

Run Individual Tests

A collection of tests is available for testing functionality of the demo adaptor. These are all contained in the subfolders of *matlabroot*\toolbox\daq\daqsdk\tests\+daq\+sdk\+tests. Each test file name begins with the letter t and has the extention .m.

To get help and information on running an individual test, use the MATLAB help command with the full package and test name. For example, to learn about the test defined in *matlabroot*\toolbox\daq\daqsdk\tests\+daq\+sdk\+tests\+workflow\tinputsinglescan.m, type:

help daq.sdk.tests.workflow.tinputsinglescan

As indicated in the display help, you can run this test with the following commands:

```
t = daq.sdk.tests.workflow.tinputsinglescan;
results = run(t);
table(results)
```

Running daq.sdk.tests.workflow.tinputsinglescan .. Done daq.sdk.tests.workflow.tinputsinglescan

Name	Passed	Failed	Incomplete	Duration
'daq.sdk.tests.workflow.tinputsinglescan/verifyInputSingleScan'	true	false	false	1.7093
'daq.sdk.tests.workflow.tinputsinglescan/verifyInputSingleScanLoop'	true	false	false	1.3631

Tip When modifying functionality in your custom adaptor, you should also modify the corresponding test. Be sure that the test runs as expected before moving on to your next modification.

Run a Test Suite

You can run all the tests in a package folder using the runtests function. For example, to run all the tests contained in daq\+sdk\+tests\+workflow, use the following commands:

```
results = runtests('daq.sdk.tests.workflow','Verbosity','Concise');
table(results)
```

Name	Passed	Failed	Incomplete
'daq.sdk.tests.workflow.tbackground/verifyAnalogInputSession'	true	false	false
'daq.sdk.tests.workflow.tbackground/verifyAnalogOutputSession'	true	false	false
'daq.sdk.tests.workflow.tbackground/verifyAnalogInputContinuous'	true	false	false
'daq.sdk.tests.workflow.tbackground/verifyAnalogOutputContinuous'	true	false	false

To run a suite of tests that includes all subpackages of a specific package, use the 'IncludeSubpackages' option in the runtests function call. The following code runs all tests below the tests package:

```
results = runtests('daq.sdk.tests','IncludeSubpackages',true,'Verbosity','Concise');
table(results)
```

Tip Run your complete modified test suite when all your individual updates are implemented and built.

See Also

Functions enableDemoAdaptorDiscovery | run | runtests

Related Examples

- "Demo Adaptor Description" on page 2-2
- "Enable the Demo Adaptor" on page 2-5
- "Create Your Adaptor from the Demo Adaptor" on page 3-2
- "Modify Demo Tests for Your Adaptor" on page 3-7

Custom Adaptor Creation

- "Create Your Adaptor from the Demo Adaptor" on page 3-2
- "Modify Demo Tests for Your Adaptor" on page 3-7
- "Errors and Exceptions" on page 3-11
- "Channel Groups" on page 3-13
- "Custom Functions" on page 3-15
- "Vendor Adaptor Templates" on page 3-17
- "Deliver Your Adaptor" on page 3-19

Create Your Adaptor from the Demo Adaptor

Use the demo adaptor as a template for creating a custom adaptor which you can build, test, and access from the toolbox. The following sections provide a sequence of steps for adaptor modification. The examples in this topic create a custom adaptor named MyAdaptor with a vendor ID of my.

In this section ...

"Edit and Build Your Adaptor" on page 3-2

"Use Your Adaptor in a Session" on page 3-4

Edit and Build Your Adaptor

This section describes the step to make a new custom adaptor based on the shipped demo adaptor. This example modifies only the names of the adaptor, vendor, and devices, without any functional changes. You build the custom adaptor in a local folder, then add the build folders to the MATLAB path. This section uses two folder locations throughout:

Location	Description
matlabroot	MATLAB installation location. This is the MATLAB used both for the building of the adaptor, and for accessing the adaptor through a data acquisition session.
C:\adaptors\daqsdk	Local file location where the new adaptor is modified and built.

- 1 Create the build area in a location of your choice. This example works with a new folder, C:\adaptors\daqsdk. Create a subfolder here called src, and within that a subfolder named daqadaptor.
- 2 Copy the folder DemoAdaptor from *matlabroot*\toolbox\daq\daqsdk\src \daqadaptor into C:\adaptors\daqsdk\src\daqadaptor.
- 3 Inside C:\adaptors\daqsdk\src\daqadaptor, rename the folder DemoAdapter to be MyAdaptor.

Navigate into MyAdaptor, and rename three of its files according to the following table:

Original Name	New Name
custom_demo.cpp	custom_my.cpp
demoadaptor.cpp	myadaptor.cpp
demoadaptor.hpp	myadaptor.hpp

- 4 With a text editor, modify each of the three new files in the previous table, replacing all occurrences of text DemoAdaptor, demoadaptor, DemoDriver, and custom_demo.cpp with MyAdaptor, myadaptor, MyDriver, and custom_my.cpp, respectively, keeping the letter capitalization style with each replacement.
- **5** Further edit the contents of myadaptor.cpp as shown in the following table:

Original Text	Updated Text
<pre>{ shortName = "MW"; fullName = "MathWorks"; driverName = "DemoAdaptor"; return DAQSuccess; }</pre>	<pre>{ shortName = "MY"; fullName = "MyAdaptor"; driverName = "MyDriver"; return DAQSuccess; }</pre>
<pre>prefix = "MWDev";</pre>	<pre>prefix = "MyDev";</pre>

With these modifications saved and in place, you are ready to build the adaptor.

6 In MATLAB, run the following utility to build the executable MEX-file for MyAdaptor:

daq.sdk.utility.mex.buildAdaptor('MyAdaptor','custom_my','C:\adaptors\daqsdk\src\daqadaptor\MyAdaptor',

The function input arguments specify the adaptor name, source code file, source file location, and where to put the built output.

Note The buildAdaptor function requires that your system be configured with Microsoft[®] Visual Studio[®] 2013 or later.

- 7 Create the folder C:\adaptors\daqsdk\+daq, and copy into it the folder +demoadaptor found at matlabroot\toolbox\daq\daqsdk\+daq\ +demoadaptor.
- 8 Navigate into C:\adaptors\daqsdk\+daq, and rename +demoadaptor to +myadaptor.

9 Navigate into C:\adaptors\daqsdk\+daq\+myadaptor, and edit these two MATLAB files in that folder:

Session.m VendorInfo.m

- In both of these files, replace all occurrences of the texts DemoAdaptor and demoadaptor with MyAdaptor and myadaptor, respectively, keeping the letter capitalization style with each replacement.
- In the VendorInfo file, use % characters to comment out the lines that hide the adaptor, between the begin and end remove indicators. The change looks like this:

```
% BEGIN REMOVE
% if daq.internal.getOptions().HideDAQSDKAdaptor
% throw(MException(message('daqsdk:HardwareInfo:VendorIsHidden', mfilename('class'))));
% end
% END REMOVE
```

• Save and close the files.

Your modified adaptor is now ready for use.

Use Your Adaptor in a Session

This example shows how to access the vendor and device represented by your modified adaptor. A data acquisition session with your adaptor allows you to add channels and get information from the device.

Start MATLAB, and use the following commands to make your adaptor available.

```
addpath 'C:\adaptors\daqsdk\bin\win64'
addpath 'C:\adaptors\daqsdk'
```

Then you can access your adaptor.

```
v = daq.getVendors
v =
Number of vendors: 2
index
        ID
              Operational
                               Comment
              -----
1
    ni
              false
                      Click here for more info
2
    mν
              true
                        MyAdaptor
```

Use the index of the vendor ID my to get more information.

```
vendor = v(2)
vendor =
Data acquisition vendor 'MyAdaptor':
           ID: 'my'
     FullName: 'MyAdaptor'
AdaptorVersion: '3.11 (R2017a)'
DriverVersion: '1.0.0'
IsOperational: true
Create a session for your device.
s = daq.createSession('my')
s =
Data acquisition session using MyAdaptor hardware:
  Will run for 1 second (1000 scans) at 1000 scans/second.
  No channels have been added.
ch1 = addAnalogOutputChannel(s, 'MyDevO', 'aoO', 'Voltage')
ch1 =
Data acquisition analog output voltage channel 'ao0' on device 'MyDevO':
TerminalConfig: SingleEnded
         Range: -10 to +10 Volts
          Name: ''
            ID: 'ao0'
        Device: [1x1 daq.sdk.DeviceInfo]
MeasurementType: 'Voltage'
s
s =
Data acquisition session using MyAdaptor hardware:
  No data queued. Will run at 1000 scans/second.
  Number of channels: 1
     index Type Device Channel MeasurementType
                                                        Range
                                                                   Name
     ---- ---
                -----
     1
           ao MyDev0 ao0
                            Voltage (SingleEnd) -10 to +10 Volts
whos
  Name
           Size
                            Bytes Class
                                                                        Attributes
  ch1
                                8 dag.sdk.AnalogOutputVoltageChannel
           1x1
                                8 dag.myadaptor.Session
  s
           1x1
           1x2
                              16 dag VendorInfo
  v
  vendor
                                8 daq.myadaptor.VendorInfo
           1x1
```

See Also

Functions buildAdaptor

Related Examples

• "Modify Demo Tests for Your Adaptor" on page 3-7

More About

• "Adaptor Creation Summary" on page 1-4

Modify Demo Tests for Your Adaptor

This topic describes how to copy demo adaptor tests and modify them for use with your own adaptor. The steps below assume you have an adaptor called MyAdaptor, as created in the example of "Create Your Adaptor from the Demo Adaptor" on page 3-2.

- 1 Copy matlabroot\toolbox\daq\daqsdk\tests to C:\adaptors\daqsdk
 \tests
- 2 In a file browser, navigate to the SDK tests package folder C:\adaptors\daqsdk \tests\+daq\+sdk.
- **3** Rename the folder +tests to +mytests.

The next steps require you to edit and save your test files. You can use the MATLAB editor, or any editor of your choice. Because the tests are MATLAB files, using the MATLAB editor is recommended for debugging purposes.

4 Navigate to C:\adaptors\daqsdk\tests\+daq\+sdk\+mytests, and open the file hardwareconfiguration.m. In MATLAB you can navigate to its folder and open the editor:

```
cd ('C:\adaptors\daqsdk\tests\+daq\+sdk\+mytests')
edit hardwareconfiguration
```

Change the vendor and device parameters in this manner, using your own names.

Original Text	Updated Text
% HardwareInfo VendorName = 'mw'; VendorFullName = 'MathWorks';	<pre>% HardwareInfo VendorName = 'my'; VendorFullName = 'MyAdaptor';</pre>
<pre>DeviceID1 = 'MWDev0'; DeviceID2 = 'MWDev1'; DeviceID3 = 'MWDev2';</pre>	DeviceID1 = 'MyDev0'; DeviceID2 = 'MyDev1'; DeviceID3 = 'MyDev2';

5 Save and close the file.

The updated vendor information now allows your tests to run on your adaptor.

6 Modify all files in C:\adaptors\daqsdk\tests\+daq\+sdk\+mytests\ +workflow\ so that all lines use mytests instead of tests. For example,

classdef tbackground < daq.sdk.mytests.workflow.BaseDAQSessionWorkflowTester</pre>

7 Restart MATLAB. Use the following commands to add your adaptor and tests to the command path.

```
addpath 'C:\adaptors\daqsdk\bin\win64'
addpath 'C:\adaptors\daqsdk'
addpath 'C:\adaptors\daqsdk\tests'
```

Run the test for single scan inputs.

```
t = daq.sdk.mytests.workflow.tinputsinglescan;
results = run(t);
table(results)
```

Running daq.sdk.mytests.workflow.tinputsinglescan

```
Done daq.sdk.mytests.workflow.tinputsinglescan
```

```
ans =
```

. .

```
2×6 table
```

Name	Passed
'daq.sdk.mytests.workflow.tinputsinglescan/verifyInputSingleScan'	true
'daq.sdk.mytests.workflow.tinputsinglescan/verifyInputSingleScanLoop'	true

8 For streaming tests, there are three files to modify in the folder C:\adaptors \daqsdk\tests\+daq\+sdk\+mytests\+development\+streaming.

Modify tstreambasic.m using your own vendor and device information, as follows:

Original Text	Updated Text	
<pre>properties(TestParameter)</pre>	<pre>properties(TestParameter)</pre>	
<pre>VendorName = {'mw'}; % Add v</pre>	<pre>VendorName = {'my'}; % Add v</pre>	endor ada
<pre>DeviceID = {'MWDev1'}; % Add</pre>	DeviceID = {'MyDev1'}; % Add	device I
end	end	

Modify both tstreamread.m and tstreamwrite.m using your own adaptor information, as follows:

Original Text	Updated Text	
<pre>properties (ClassSetupParameter) adaptorName = { 'DemoAdaptor' }</pre>	<pre>properties (ClassSetupParameter) adaptorName = {'MyAdaptor'}</pre>	
end	end	
<pre>adaptorPath = fullfile(matlabroot,'t</pre>	adaptorPath = 'c:\adaptors\daqsdk\bi	rn(\\waiumc6h4'');

9 You can now run any of the streaming tests on your adaptor. For example, restart MATLAB and enter the following code:

```
addpath 'C:\adaptors\daqsdk\bin\win64'
addpath 'C:\adaptors\daqsdk'
addpath 'C:\adaptors\daqsdk\tests'
t = daq.sdk.mytests.development.streaming.tstreamread;
results = run(t);
table(results)
```

Running daq.sdk.mytests.development.streaming.tstreamread Done daq.sdk.mytests.development.streaming.tstreamread

ans =

8×6 table

Name

'daq.sdk.mytests.development.streaming.tstreamread[adaptorName=MyAdaptor]/Verif 'daq.sdk.mytests.development.streaming.tstreamread[adaptorName=MyAdaptor]/Verif 'daq.sdk.mytests.development.streaming.tstreamread[adaptorName=MyAdaptor]/Verif 'daq.sdk.mytests.development.streaming.tstreamread[adaptorName=MyAdaptor]/Verif 'daq.sdk.mytests.development.streaming.tstreamread[adaptorName=MyAdaptor]/Verif 'daq.sdk.mytests.development.streaming.tstreamread[adaptorName=MyAdaptor]/Verif 'daq.sdk.mytests.development.streaming.tstreamread[adaptorName=MyAdaptor]/Verif 'daq.sdk.mytests.development.streaming.tstreamread[adaptorName=MyAdaptor]/Verif 'daq.sdk.mytests.development.streaming.tstreamread[adaptorName=MyAdaptor]/Verif

Further Suggestions

Run Test Suites

You can run the full suite of tests for your adaptor by specifying the package folder to use all the tests contained in it.

```
results = runtests('daq.sdk.mytests','IncludeSubpackages',true,'Verbosity','Concise');
table(results)
```

Modify Functionality Tests

- As you write your adaptor, you must modify the test files to correspond to the functionality implemented for your device. In deciding the sequence in which you implement and test functionality, consider "Session Workflows with the Demo Adaptor" on page 2-6.
- In addition to adaptor name, you must modify where applicable the vendor name, device driver name, device names, vendor ID, etc.

See Also

Functions run | runtests

Related Examples

- "Test the Demo Adaptor" on page 2-12
- "Create Your Adaptor from the Demo Adaptor" on page 3-2

Errors and Exceptions

Nonstreaming

To indicate that a standard SDK function has resulted in an error for an expected reason, return the appropriate error code (as opposed to DAQSuccess), as provided in include/ daqsdktypes.h.

To indicate that a custom SDK function has resulted in an error, throw a DAQDiagnostic (see daqinterfaces.hpp) containing a custom error code and a diagnosis message string.

To indicate that a standard SDK function has resulted in an error for a reason that is specific to the function of the custom adaptor (vendor-specific error), throw a DAQDiagnostic.

You can define a custom error code as a negative value less than daqsdk::DAQErr_ReservedRangeEnd (see daqsdktypes.h). while a custom warning code can be defined as a positive value greater than daqsdk::DAQWrn_ReservedRangeEnd.

Streaming

To indicate that an error has occurred during the configuration of the stream (configureStream), return a custom error code. You should also implement the DAQStreamAnalog::getDiagnosticFromStatus, which when given a custom error code, returns a string describing the error condition.

To indicate that an error has occurred during streaming (that is, after the stream has started but before it is done or has stopped), return a custom error code or throw an exception.

Related Examples

• "Create Your Adaptor from the Demo Adaptor" on page 3-2

More About

• "Adaptor API Reference" on page 4-2

- "Streaming API Reference" on page 4-21
- "Sequence for Errors and Exceptions" on page 5-21

Channel Groups

In this section ...

"Channel Group Description" on page 3-13

"Channel Group Restrictions" on page 3-14

Channel Group Description


A simplified view of a DAQ device is that actual devices provide channels with common functions, logically grouped into subsystems. For example, all channels that provide analog input data may be thought of as belonging to an analog input subsystem.

Another view of this DAQ device is as a provider and consumer of data. The largest unit of data that can be acquired or generated simultaneously, by one or more channels, is a scan. The logical grouping that acquires or consumes one or more scans of data is a channel group. The definition of the channel group is usually constrained by the driver and hardware. For example, when all channels belonging to a single analog input subsystem also share a single clock.

A *channel group* is an aggregation of channels, usually of the same subsystem, which operate together. For example, all the analog output subsystem channels on a device must be configured, reserved, and act together to generate data as a single scan.

You should define channel groups in a way that reflects the driver constraints, and provide a means for identifying all channels belonging to the group for acquiring and generating scans. Typically, channel groups provide functions to stream data to or from a device buffer.

Each channel on the device has a unique address, defined by device, subsystem, and channel ID. Each channel must be assigned to one channel group. The following diagram illustrates one possible channel group arrangement. In this scenario, analog output (AO) channels 0 and 1 might serve a different purpose than AO channels 2 and 3; while analog input (AI) channels 0–3 are used all together.

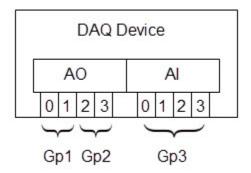

For reference information on the functions used in configuring channel groups, see "Hardware Management" on page 4-7.

Channel Group Restrictions

- All channels in a group operate together. This allows synchronized streaming to the extent supported by the hardware.
- A channel cannot belong to more than one group.
- · All channels in a group are requested, reserved, and released together.
- By default, a channel group—and therefore all its channels—can be accessed by only one data acquisition session at a time. You cannot add channels to a session if any other channels in their groups are already added to a different session. If your driver allows a group to be accessed by different sessions, you can control this behavior using the isRegistrationReservationImpl function.

More About

- "Vendor Adaptor Templates" on page 3-17
- "Adaptor API Reference" on page 4-2
- "Streaming API Reference" on page 4-21
- "Streaming Input and Output" on page 2-10
- "Streaming Sequence Diagrams" on page 5-4

Custom Functions

Custom functionality provided by your adaptor that is not part of the standard sessionbased interface can be exposed to MATLAB via the DAQ SDK custom interface. For example, your device might provide an on-board power supply.

Note This topic assumes experience writing MEX-files.

To add a custom function, first review the custom functions available in the demo adaptor in the folder *matlabroot*\matlab\toolbox\daq\daqsdk\+daq\ +demoadaptor\+custom. The installed files in this folder are:

testHasInputsHasOutputs.m
testHasInputsNoOutputs.m
testNoInputsNoOutputs.m
testThrowCustomExceptions.m

Use these steps to create your own custom function:

- 1 Add a function to the MyDriver class (MyDriver::customFunction).
- 2 Add a function to the MyAdaptor class (MyAdaptor::customFunction) that calls MyDriver::customFunction with the designated:
 - Inputs
 - Outputs
 - Custom error code
- **3** Update custom_my.cpp to:
 - Define a function to call (dispatch) the custom adaptor function MyAdaptor::customFunction.
 - Update the customizeMap function to add:

functionMap["myCustomFunction"] = customFunction;

where customFunction is the name of the MEX-function that calls MyAdaptor::customFunction, and myCustomFunction is the name of function in MATLAB.

- 4 Define your custom MATLAB function myCustomFunction.m in the +daq\ +myadaptor\+custom subpackage for your adaptor.
 - Choose the appropriate template from \+daq\+demoadaptor\+custom\ to copy and rename.
 - Has Inputs, Has Outputs
 - Has Inputs, No Outputs
 - No Inputs, No Outputs
 - Rename the file to perform the desired function, for example, myCustomFunction.m.
 - Edit myCustomFunction.m to
 - Update: functionName = 'myCustomFunction';
 - Provide the inputs to the function as a structure.

For functionality that is not part of the standard session interface, contact MathWorks[®] technical support at https://www.mathworks.com/support/contact_us to let us know what functionality you need.

More About

- "C/C++ Matrix Library API" (MATLAB)
- "Vendor Adaptor Templates" on page 3-17
- "Errors and Exceptions" on page 3-11
- "Deliver Your Adaptor" on page 3-19

Vendor Adaptor Templates

The Data Acquisition Toolbox SDK is installed with a set of source file stubs for an adaptor called VendorAdaptor. The source files are installed in the folder:

matlabroot\toolbox\daq\daqsdk\src\daqadaptor\VendorAdaptor

The basic components and structure of the source files for this adaptor are the same as those in the DemoAdaptor. If you need to create an adaptor from scratch, it is recommended that you use a copy of the VendorAdaptor source files. The following table indicates the purpose of each source file.

File	Description
vendoradaptor.hpp, vendoradaptor.cpp	Wraps device driver code in methods that allow you to configure, discover, and enumerate the hardware in MATLAB.
daqstream_analog.hpp, daqstream_analog.cpp	Implements a DAQStream object for an analog subsystem, that allows you to stream data to and from the hardware.
custom_vendor.cpp	Dispatches custom calls from MATLAB to the adaptor. At a minimum this must contain a customizeMap function.

Typical Workflow to Create Adaptor

To create an adaptor from the set of template files in the folder VendorAdaptor, use the following steps. Assume that you want to name the adaptor MyAdaptor.

- 1 Create a copy of the entire folder, and name it MyAdaptor.
- 2 Working in the new folder called MyAdaptor, change the names of the files:

Original Name	New Name
vendoradaptor.hpp	myadaptor.hpp
vendoradaptor.cpp	myadaptor.cpp
custom_vendor.cpp	custom_my.cpp

3 Update the content of the files so that the new names are used for references to other files, the adaptor, devices, and vendor.

- **4** Update the functions to use your driver code. For more information, see "Adaptor API Reference" on page 4-2.
- **5** Build the adaptor with the buildAdaptor function.

Related Examples

• "Create Your Adaptor from the Demo Adaptor" on page 3-2

More About

- "Adaptor Creation Summary" on page 1-4
- "Custom Functions" on page 3-15

Deliver Your Adaptor

When you have a custom adaptor for delivery, you can create a toolbox to deliver the adaptor MEX-file along with your device driver and any other support files, such as documentation, data files, examples, and so on.

You should document any adaptor behavior that involves:

- · Differences from standard data acquisition session behavior
- Custom functions

For information on creating and distributing a toolbox, see "Toolbox Distribution" (MATLAB).

For creating and delivering documentation to support your toolbox, see "Display Custom Documentation" (MATLAB).

For your custom examples, see "Display Custom Examples" (MATLAB).

API Reference

- "Adaptor API Reference" on page 4-2
- "Streaming API Reference" on page 4-21

Adaptor API Reference

This topic provides an overview of each function included in the demo adaptor source file, demoadaptor.cpp, grouped in the following categories. The vendoradaptor.cpp template includes similar functions.

In this section...

"Lifetime" on page 4-2 "Enumeration" on page 4-3 "Hardware Management" on page 4-7 "Vendor and Device Discovery" on page 4-8 "Subsystem Discovery" on page 4-10 "Configuration" on page 4-15 "Reservation" on page 4-18 "SingleScans" on page 4-19

Lifetime

Lifetime functions include those that involve the loading and unloading of the driver interface.

initImpl

Syntax	DAQStatus DemoDriver::initImpl()
Purpose	Initialize and load implementation of daqsdk::IDriver interface
Inputs	None
Output	None
Return status	DAQErr_Driver_Init on failure.
	DAQSuccess on success.

termImpl

Syntax	DAQStatus DemoDriver::termImpl()
Purpose	Terminate and unload implementation of daqsdk::IDriver interface

Inputs	None
Output	None
Return status	DAQErr_Driver_Term on failure.
	DAQSuccess on success.

Enumeration

Enumeration functions involve the recognition of devices, subsystems, and channels.

enumerateDevicesImpl

Syntax	<pre>DemoDriver::enumerateDevicesImpl(Index &deviceCount) const</pre>
Purpose	Enumerate devices available via vendor driver
Inputs	None
Output	Number of the devices enumerated
Return status	DAQErr_Driver_EnumerateDevices on failure.
	DAQSuccess on success.

commitDevicesImpl

Syntax	<pre>DemoDriver::commitDevicesImpl(Index deviceCount)</pre>
Purpose	Inform the driver that the enumerated devices are to be committed, in enumerated order, for use by the adaptor
Inputs	Number of devices enumerated
Output	None
Return status	DAQErr_Driver_CommitDevices on failure.
	DAQSuccess on success.

enumerateSubsystemsImpl

Syntax	<pre>DemoDriver::enumerateSubsystemsImpl(Index deviceIndex,</pre>	
	Index &subsystemCount) const	

Purpose	Enumerate the subsystems available via for a given device
Inputs	Index of the given device
Output	Number of the subsystems enumerated
Return status	DAQErr_Driver_EnumerateSubsystems on failure.
	DAQSuccess on success.

commitSubsystemsImpl

Syntax	<pre>DemoDriver::commitSubsystemsImpl(Index deviceIndex, Index subsystemCount)</pre>
Purpose	Inform the driver that the enumerated subsystems, for a given device, are to be committed, in enumerated order, for use by the adaptor
Inputs	Index of the given device, number of the subsystems enumerated
Output	None
Return status	DAQErr_Driver_CommitSubsystems on failure.
	DAQSuccess on success.

enumerateChannelsImpl

Syntax	DemoDriver::enumerateChannelsImpl(Index deviceIndex, Index subsystemIndex, Index &channelCount) const
Purpose	Enumerate the channels available via for a given subsystem and device
Inputs	Index of the given device, the index of the given subsystem
Output	Number of the channels enumerated
Return status	DAQErr_Driver_EnumerateChannels on failure.
	DAQSuccess on success.

commitChannelsImpl

Syntax	<pre>DemoDriver::commitChannelsImpl(Index deviceIndex, Index</pre>
	subsystemIndex, Index channelCount)

Purpose	Inform the driver that the enumerated channels, for a given device and subsystem, are to be committed, in enumerated order, for use by the Adaptor
Inputs	Index of the given device, the index of the given subsystem, the number of channels enumerated
Output	None
Return status	DAQErr_Driver_CommitChannels on failure.
	DAQSuccess on success.

getChannelGroupIndexImpl

DemoDriver::getChannelGroupIndexImpl(Index deviceIndex, Index subsystemIndex, Index channelIndex, Index &channelGroupIndex) const
Return the channel group index corresponding to a specified channel
Index of the specified device, index of the specified subsystem, index of the specified channel
Channel group index
DAQErr_Driver_GetChannelGroupIndex on failure. DAQSuccess on success.

determineOrderOfChannelAdditionImpl

	DemoDriver::getOrderOfChannelAdditionImpl(daqsdk::OrderOf &orderOfChannelsInGroup)	Channels
-	Return an enumeration representing the order in which channels indices are sorted, by the driver, in channel groups	
Inputs	None	
Output	Order of channels within channel groups	
Return status	DAQErr_Driver_OrderOfChannelAddition on failure.	
	DAQSuccess on success.	

Channel groups contain a list of channels ordered first by device, then by subsystem, and finally by channel. The group must acquire data from requested channels either in the

listed order ("Sorted") or in the order requested ("InOrderOfAddition"). For example, if the group contains four channels and a user requests channels 4, 2, and 1, they should expect data from the channel group either in the order 4, 2, 1 (the order in which the channels were added) or in the order 1, 2, 4 (sorted). See "Channel Groups" on page 3-13.

releaseDevicesImpl

Syntax	<pre>DemoDriver::releaseDevicesImpl(Index deviceIndex)</pre>
Purpose	Release the resources committed by the driver for a specified device
Inputs	Index of the device resources to release
Output	None
Return status	DAQErr_Driver_ReleaseDevices on failure.
	DAQSuccess on success.

releaseSubsystemsImpl

Syntax	<pre>DemoDriver::releaseSubsystemsImpl(Index deviceIndex, Index subsystemIndex)</pre>
Purpose	Release the resources committed by the driver for a specified subsystem and device
Inputs	Index of the device resources to release, index of the subsystem resources to release
Output	None
Return status	DAQErr_Driver_ReleaseSubsystems on failure. DAQSuccess on success.

releaseChannelsImpl

Syntax	<pre>DemoDriver::releaseChannelsImpl(Index deviceIndex, Index subsystemIndex, Index channelIndex)</pre>
Purpose	Release the resources committed by the driver for a specified channel of a subsystem of a device

-	Index of the device resources to release, index of the subsystem resources to release, index of the channel resources to release
Output	None
Return status	DAQErr_Driver_ReleaseChannels on failure.
	DAQSuccess on success.

Hardware Management

Hardware management functions control the configuration of channel groups.

addChannelImpl

Syntax	DemoDriver::addChannelImpl(Index deviceIndex, Index subsystemIndex, Index channelIndex)
Purpose	Register the specified channel with its channel group
Inputs	Index of the device, index of the subsystem for given device, index of the channel for the given subsystem
Output	None
Return status	DAQErr_Driver_AddChannel on failure.
	DAQSuccess on success.

removeChannelImpl

Syntax	DemoDriver::removeChannelImpl(Index deviceIndex, Index subsystemIndex, Index channelIndex)
Purpose	Unregister the specified channel from its channel group
Inputs	Index of the device, index of the subsystem for given device, index of the channel for the given subsystem
Output	None
Return status	DAQErr_Driver_RemoveChannel on failure.
	DAQSuccess on success.

reserveChannelGroupImpl

Syntax	<pre>DemoDriver::reserveChannelGroupImpl(ChannelGroupIndex groupIndex)</pre>
Purpose	Reserve the specified channel group and all its resources
Inputs	Index of the channel group
Output	None
Return status	DAQErr_Driver_ReserveChannelGroup on failure.
	DAQSuccess on success.

unreserveChannelGroupImpl

Syntax	<pre>DemoDriver::unreserveChannelGroupImpl(ChannelGroupIndex groupIndex)</pre>
Purpose	Unreserve/release the specified channel group and all its resources
Inputs	Index of the channel group
Output	None
Return status	DAQErr_Driver_UnreserveChannelGroup on failure.
	DAQSuccess on success.

Vendor and Device Discovery

These functions retrieve information about vender and device.

getDriverVersionImpl

Syntax	DemoDriver::getDriverVersionImpl(uint32_T &major, uint32_T &minor, uint32_T &patch) const
Purpose	Return driver version number
Inputs	None
Output	Major version number, minor version number, patch version number

Return status	DAQErr_Driver_GetDriverVersion on failure.
	DAQSuccess on success.

getVendorInfoImpl

Syntax	<pre>DemoDriver::getVendorInfoImpl(std::string &shortName, std::string &fullName, std::string &driverName) const</pre>
Purpose	Return relevant vendor information (name and driver-name)
Inputs	None
Output	Vendor shortname (typically used as a vendor ID), vendor fullname, driver name (including full path)
Return status	DAQErr_Driver_GetVendorInfo on failure.
	DAQSuccess on success.

getDeviceInfoImpl

Syntax	<pre>DemoDriver::getDeviceInfoImpl(Index deviceIndex, std::string &model, std::string &prefix, std::string &id, std::string &serialNumber, bool &isRecognizedDevice) const</pre>
Purpose	Return relevant device information
Inputs	Index of the device
Output	Device model, device prefix (e.g., 'Dev', 'Audio', etc.), device ID, device serial number, indication of whether the driver recognizes and supports the device
Return status	DAQErr_Driver_GetDeviceInfo on failure.
	DAQSuccess on success.

getFirmwareVersionImpl

Syntax	<pre>DemoDriver::getFirmwareVersionImpl(Index deviceIndex,</pre>
	uint32_T &major, uint32_T &minor, uint32_T &patch) const
Purpose	Return firmware version number

Inputs	None
Output	Major version number, minor version number, patch version number
Return status	DAQErr_Driver_GetFirmwareVersion on failure.
	DAQSuccess on success.

Subsystem Discovery

These functions retrieve information about the subsystem.

getSubsystemsOfTypeImpl

Syntax	<pre>DemoDriver::getSubsystemsOfTypeImpl(Index deviceIndex, IndexList &subsystemIndices, daqsdk::Subsystem subsystemType, daqsdk::TransferDirection transferDirection) const</pre>
Purpose	Return subsystems of a given type (Analog, Digital, etc.) and direction (Input, Output)
Inputs	Index of the device, subsystem type, transfer direction
Output	List of subsystem indices with the given type/direction or empty if no matches are found
Return status	request (but not when no subsystems are found).
	DAQSuccess on success.

getMeasurementTypesImpl

Syntax	<pre>DemoDriver::getMeasurementTypesImpl(Index deviceIndex, Index subsystemIndex, std::vector<daqdatatypes::measurementtype> &measurementTypes) const</daqdatatypes::measurementtype></pre>
Purpose	Return the measurement types supported by a specified subsystem and device
Inputs	Index of the device, index of the subsystem
Output	List of measurement types supported by the specified subsystem

	DAQErr_Driver_GetMeasurementTypes on failure to execute the request (but not when no subsystems are found).
	DAQSuccess on success.

getDefaultMeasurementTypeImpl

Syntax	<pre>DemoDriver::getDefaultMeasurementTypeImpl(Index deviceIndex, Index subsystemIndex, daqdatatypes::MeasurementType &defaultMeasurementType) const</pre>
Purpose	Return the default measurement type supported by a specified subsystem and device
Inputs	Index of the device, index of the subsystem
Output	Default measurement types supported by the specified subsystem
Return status	
	DAQSuccess on success.

getRateLimitImpl

Syntax	<pre>DemoDriver::getRateLimitImpl(Index deviceIndex, Index subsystemIndex, daqdatatypes::RateLimit &rateLimit) const</pre>
Purpose	Return the rate limits supported by a specified subsystem and device
Inputs	Index of the device, index of the subsystem
Output	Rate limits supported by the specified subsystem
Return status	DAQErr_Driver_GetRateLimit on failure.
	DAQSuccess on success.

getResolutionImpl

Syntax	<pre>DemoDriver::getResolutionImpl(Index deviceIndex, Index subsystemIndex, uint8_T &resolution) const</pre>
Purpose	Return the measurement resolution supported by a specified subsystem and device

Inputs	Index of the device, index of the subsystem
Output	Measurement resolution supported by the specified subsystem
Return status	DAQErr_Driver_GetResolution on failure.
	DAQSuccess on success.

getTerminalConfigsImpl

Syntax	<pre>DemoDriver::getTerminalConfigsImpl(Index deviceIndex, Index subsystemIndex, std::vector<daqdatatypes::terminalconfiguration> &terminalConfigurations) const</daqdatatypes::terminalconfiguration></pre>
Purpose	Return the terminal configurations supported by a specified subsystem and device for each channel
Inputs	Index of the device, index of the subsystem
Output	Terminal configurations supported by the specified subsystem
Return status	DAQErr_Driver_GetTerminalConfigs on failure.
	DAQSuccess on success.

getRangesAvailableForTerminalConfigImpl

Syntax	<pre>DemoDriver::getRangesAvailableForTerminalConfigImpl(Index deviceIndex, Index subsystemIndex, daqdatatypes::TerminalConfiguration terminalConfig, std::vector<daqdatatypes::range> &ranges) const</daqdatatypes::range></pre>
Purpose	Return ranges supported by specified terminal configuration for subsystem and device
Inputs	Index of the device, index of the subsystem, terminal configuration type
Output	Ranges supported by the specified terminal configuration for a given subsystem
Return status	DAQErr_Driver_GetRangesAvailableForTerminalConfig on failure.
	DAQSuccess on success.

getDefaultTerminalConfigsImpl

Syntax	<pre>DemoDriver::getDefaultTerminalConfigsImpl(Index deviceIndex, Index subsystemIndex, std::vector<daqdatatypes::terminalconfiguration> &defaultTerminalConfigs) const</daqdatatypes::terminalconfiguration></pre>
Purpose	Return the default terminal configuration types supported by a specified subsystem and device
Inputs	Index of the device, index of the subsystem
Output	Default terminal configuration types supported by the specified subsystem
Return status	DAQErr_Driver_GetDefaultTerminalConfigs on failure.
	DAQSuccess on success.

is On Demand Operation Supported Impl

Syntax	<pre>DemoDriver::isOnDemandOperationSupportedImpl(Index deviceIndex, Index subsystemIndex, bool &isSupported) const</pre>
Purpose	Indicate whether on-demand operations are supported by a specified subsystem and device
Inputs	Index of the device, index of the subsystem
Output	Whether on-demand operations are supported by the specified subsystem
Return status	DAQErr_Driver_IsOnDemandOperationSupported on failure.
	DAQSuccess on success.

getCouplingsImpl

Syntax	<pre>DemoDriver::getCouplingsImpl(Index deviceIndex, Index subsystemIndex, std::vector<daqdatatypes::coupling> &couplings) const</daqdatatypes::coupling></pre>
Purpose	Return the couplings supported by a specified subsystem and device
Inputs	Index of the device, index of the subsystem

Output	Couplings supported by the specified subsystem
Return status	DAQErr_Driver_GetCouplings on failure.
	DAQSuccess on success.

getDefaultCouplingImpl

Syntax	DemoDriver::getDefaultCouplingImpl(Index deviceIndex, Index subsystemIndex, daqdatatypes::Coupling &defaultCoupling) const
Purpose	Return the default coupling supported by a specified subsystem and device
Inputs	Index of the device, index of the subsystem
Output	Default coupling supported by the specified subsystem
Return status	DAQErr_Driver_GetDefaultCoupling on failure.
	DAQSuccess on success.

getSampleTypesImpl

Syntax	<pre>DemoDriver::getSampleTypesImpl(Index deviceIndex, Index subsystemIndex, std::vector<daqdatatypes::sampletype> &sampleTypes) const</daqdatatypes::sampletype></pre>
Purpose	Return the sample types supported by a specified subsystem and device
Inputs	Index of the device, index of the subsystem
Output	Sample types supported by the specified subsystem
Return status	DAQErr_Driver_GetSampleTypes on failure.
	DAQSuccess on success.

getDefaultSamplingTypeImpl

Syntax	DemoDriver::getDefaultSamplingTypeImpl(Index
	deviceIndex, Index subsystemIndex,
	daqdatatypes::SampleType &defaultSampleType)

Purpose	Return the default sample type supported by a specified subsystem and device
Inputs	Index of the device, index of the subsystem
Output	Default sample type supported by the specified subsystem
Return status	DAQErr_Driver_GetDefaultSamplingType on failure.
	DAQSuccess on success.

getChannelNamesImpl

Syntax	<pre>DemoDriver::getChannelNamesImpl(Index deviceIndex, Index subsystemIndex, std::vector<std::string> &channelNames) const</std::string></pre>
Purpose	Return the channel names supported by a specified subsystem and device
Inputs	Index of the device, index of the subsystem
Output	Channel names supported by the specified subsystem
Return status	DAQErr_Driver_GetChannelNames on failure.
	DAQSuccess on success.

Configuration

Configuration functions control rates, ranges, and coupling.

getRateImpl

Syntax	<pre>DemoDriver::getRateImpl(ChannelGroupIndex groupIndex, daqsdk::float64 &rate) const</pre>
Purpose	Return the rate supported by a specified channel group in its current configuration
Inputs	Index of the group
Output	Rate supported by the specified channel group
Return status	DAQErr_Driver_GetRate on failure.
	DAQSuccess on success.

setRateImpl

Syntax	<pre>DemoDriver::setRateImpl(ChannelGroupIndex groupIndex, daqsdk::float64 rate)</pre>
Purpose	Set the rate for a specified channel group in its current configuration
Inputs	Index of the group, rate
Output	None
Return status	DAQErr_Driver_SetRate on failure.
	DAQSuccess on success.

getChannelCouplingImpl

Syntax	DemoDriver::getChannelCouplingImpl(Index deviceIndex, Index subsystemIndex, Index channelIndex, daqdatatypes::Coupling &coupling) const
Purpose	Return the channel coupling of a specified channel for a given subsystem and device
Inputs	Index of the device, index of the subsystem, index of the channel
Output	Coupling supported by the specified channel
Return status	DAQErr_Driver_GetChannelCoupling on failure.
	DAQSuccess on success.

setChannelCouplingImpl

Syntax	<pre>DemoDriver::setChannelCouplingImpl(Index deviceIndex, Index subsystemIndex, Index channelIndex, std::string coupling)</pre>
Purpose	Set the channel coupling of a specified channel for a given subsystem and device
Inputs	Index of the device, index of the subsystem, index of the channel
Output	None
Return status	DAQErr_Driver_SetChannelCoupling on failure.
	DAQSuccess on success.

getChannelTerminalConfigImpl

Syntax	<pre>DemoDriver::getChannelTerminalConfigImpl(Index deviceIndex, Index subsystemIndex, Index channelIndex, daqdatatypes::TerminalConfiguration &terminalConfig) const</pre>
Purpose	Return the terminal configuration of a specified channel for a given subsystem and device
Inputs	Index of the device, index of the subsystem, index of the channel
Output	Terminal configuration supported by the specified channel
Return status	DAQErr_Driver_GetChannelTerminalConfig on failure.
	DAQSuccess on success.

setChannelTerminalConfigImpl

Syntax	<pre>DemoDriver::setChannelTerminalConfigImpl(Index deviceIndex, Index subsystemIndex, Index channelIndex, std::string terminalConfig)</pre>
Purpose	Set the terminal configuration of a specified channel for a given subsystem and device
Inputs	Index of the device, index of the subsystem, index of the channel
Output	None
Return status	DAQErr_Driver_SetChannelTerminalConfig on failure.
	DAQSuccess on success.

getChannelRangeImpl

Syntax	DemoDriver::getChannelRangeImpl(Index deviceIndex, Index subsystemIndex, Index channelIndex, daqdatatypes::Range ⦥) const
Purpose	Return the range of a specified channel for a given subsystem and device
Inputs	Index of the device, index of the subsystem, index of the channel
Output	Range supported by the specified channel

Return status	DAQErr_Driver_GetChannelRange on failure.
	DAQSuccess on success.

setChannelRangeImpl

Syntax	<pre>DemoDriver::setChannelRangeImpl(Index deviceIndex, Index subsystemIndex, Index channelIndex, daqdatatypes::Range range)</pre>
Purpose	Set the range of a specified channel for a given subsystem and device
Inputs	Index of the device, index of the subsystem, index of the channel
Output	None
Return status	DAQErr_Driver_SetChannelRange on failure.
	DAQSuccess on success.

Reservation

Reservation functions query device and channel availability.

is Device Available Impl

Syntax	<pre>DemoDriver::isDeviceAvailableImpl(Index deviceIndex, bool &isDeviceAvailable) const</pre>
Purpose	Return whether or not a specified device is still available, connected, committed, and enumerated
Inputs	Index of the device
Output	Return whether or not the device is available as previously committed and enumerated
Return status	DAQErr_Driver_IsDeviceAvailable on failure.
	DAQSuccess on success.

isRegistrationReservationImpl

Syntax	DemoDriver::isRegistrationReservationImpl(Index
	deviceIndex, bool &isReservation) const

Purpose	Return whether or not registering a channel in a channel group reserves the channel group
Inputs	Index of the channel group
Output	Return whether or not "registration is reservation." (See "Channel Groups" on page 3-13.)
Return status	DAQErr_Driver_IsRegistrationReservation on failure.
	DAQSuccess on success.

SingleScans

Single scan functions acquire or generate a static scan of data.

inputSingleScan

Syntax	DemoDriver::inputSingleScanImpl(ChannelGroupIndex groupIndex, DataScan &data) const
Purpose	Acquire a scan of data for all channels registered with a channel group
Inputs	None
Output	Acquired scan of data
Return status	DAQErr_Driver_InputSingleScan on failure.
	DAQSuccess on success.

OutputSingleScan

Syntax	DemoDriver::outputSingleScanImpl(ChannelGroupIndex groupIndex, DataScan &&outputData) const
Purpose	Generate a scan of data for all channels registered with a channel group
Inputs	Data to generate output
Output	None
Return status	DAQErr_Driver_OutputSingleScan on failure.
	DAQSuccess on success.

Related Examples

• "Create Your Adaptor from the Demo Adaptor" on page 3-2

More About

- "Errors and Exceptions" on page 3-11
- "Channel Groups" on page 3-13
- "Streaming API Reference" on page 4-21
- "Custom Functions" on page 3-15

Streaming API Reference

In this section...

"Initialization and Configuration" on page 4-21 "Start and Stop" on page 4-22 "Data Availability" on page 4-23 "Transfer Data" on page 4-24

Streaming is used in the generation or acquisition of clocked data to allow asynchronous operation that does not block MATLAB. Stream channels accommodate the flow of data separately from the session-dispatcher route. This also allows data sets that might exceed the size of the memory on the device.

The following functions are defined in the streaming source file daqstream_analog.cpp.

Initialization and Configuration

initialize

Syntax	<pre>int64_T DAQStreamAnalog::initialize()</pre>
Purpose	Initialize channel group after constructor.
Inputs	None
Output	None

terminate

Syntax	<pre>int64_T DAQStreamAnalog::terminate()</pre>
Purpose	Terminate DAQ stream prior to its destruction.
Inputs	None
Output	None

configureStream

Syntax	<pre>int64_T DAQStreamAnalog::configureStream()</pre>
--------	---

Purpose	Configure group of channels on the DAQ device driver for streaming operation.
Inputs	None
Output	None

unconfigureStream

Syntax	<pre>int64_T DAQStreamAnalog::unconfigureStream()</pre>
Purpose	Unconfigure groups of channels when DAQStream channel is closed from MATLAB.
Inputs	None
Output	None

registerCallbacks

Syntax	<pre>int64_T DAQStreamAnalog::registerCallbacks()</pre>
Purpose	Register any callback handlers required by DAQ device driver following <code>configureStream</code> .
Inputs	None
Output	None

unregisterCallbacks

Syntax	<pre>int64_T DAQStreamAnalog::unregisterCallbacks()</pre>
Purpose	Unregister immediately prior to unconfigureStream any callback handlers registered with DAQ device driver.
Inputs	None
Output	None

Start and Stop

prestart

Syntax	<pre>int64_T DAQStreamAnalog::prestart()</pre>
--------	--

	Called on a per-run basis prior to the streaming operation start, typically to reset scan counters.
Inputs	None
Output	None

start

Syntax	<pre>int64_T DAQStreamAnalog::start()</pre>
Purpose	Start the streaming operation for the given channelGroupHandle.
Inputs	None
Output	None

stop

Syntax	<pre>int64_T DAQStreamAnalog::stop()</pre>
Purpose	Stop the streaming operation for the given channelGroupHandle.
Inputs	None
Output	None

Data Availability

getNumInputScansAvailable

Syntax	int64_T DAQStreamAnalog::getNumInputScansAvailable(uint64_T& numScansAcquired)
Purpose	Query the number of input scans available to be read by a read or readWrite call.
Inputs	None
Output	Number of scans.

get Num Scans Output By Hardware

Syntax	int64_T
	DAQStreamAnalog::getNumScansOutputByHardware(uint64_T&
	numScansGenerated)

Purpose	Query the number of scans output by the hardware by a write or readWrite call.
Inputs	None
Output	Number of scans.

getOutputBufferSize

Syntax	<pre>int64_T DAQStreamAnalog::getOutputBufferSize(uint64_T& outputBufferSize)</pre>
Purpose	Query the DAQ device output buffer size in number of scans.
Inputs	None
Output	Buffer size in scans.

flushOutputBuffer

Syntax	<pre>int64_T DAQStreamAnalog::flushOutputBuffer()</pre>
Purpose	Empty the output buffer.
Inputs	None
Output	None

isDeviceDone

Syntax	<pre>int64_T DAQStreamAnalog::isDeviceDone(bool& isDone)</pre>
Purpose	Poll the vendor driver immediately following a call to stop.
Inputs	None
Output	True if device is done streaming.

Transfer Data

read

Syntax	int64_T DAQStreamAnalog::read(float64 * const pReadBuffer, uint64_T numReadScans)
Purpose	Read acquired data from the DAQ device into the read buffer.

Inputs	<pre>pReadBuffer: buffer used by stream to store input data acquired from the device. numReadScans: number of scans to copy into the provided buffer.</pre>
Output	None
Notes	The stream is responsible for the lifetime of the buffer.

write

Syntax	<pre>int64_T DAQStreamAnalog::write(float64 const * const pWriteBuffer, uint64_T numWriteScans)</pre>
Purpose	Write data from the buffer to the device for output generation.
Inputs	<pre>pWriteBuffer: buffer used by stream to store output data to be generated by the device. numWriteScans: number of valid scans to copy from the provided buffer.</pre>
Output	None
Notes	The stream is responsible for the lifetime of the buffer.

readWrite

Syntax	<pre>int64_T DAQStreamAnalog::readWrite(float64* const pReadBuffer, uint64_T numReadScans, float64 const * const pWriteBuffer, uint64_T numWriteScans)</pre>
Purpose	Simultaneously read and write data between the buffers and a DAQ device duplex channel.
Inputs	<pre>pReadBuffer: buffer used by stream to store input data acquired from the device. numReadScans: number of scans to copy into the provided buffer. pWriteBuffer: buffer used by stream to store output data to be generated by the device. numWriteScans: number of valid scans to copy from the provided buffer.</pre>
Output	None
Notes	The stream is responsible for the lifetime of the buffer.

Related Examples

• "Create Your Adaptor from the Demo Adaptor" on page 3-2

More About


- "Errors and Exceptions" on page 3-11
- "Channel Groups" on page 3-13
- "Adaptor API Reference" on page 4-2
- "Custom Functions" on page 3-15

State and Sequence Diagrams

- "State Machine Diagram" on page 5-2
- "Streaming Sequence Diagrams" on page 5-4
- "Foreground Streaming Sequences" on page 5-5
- "Background Streaming Sequences" on page 5-12
- "Sequence for Errors and Exceptions" on page 5-21

State Machine Diagram

This composite diagram shows the state machine for input, output, and duplex channels.

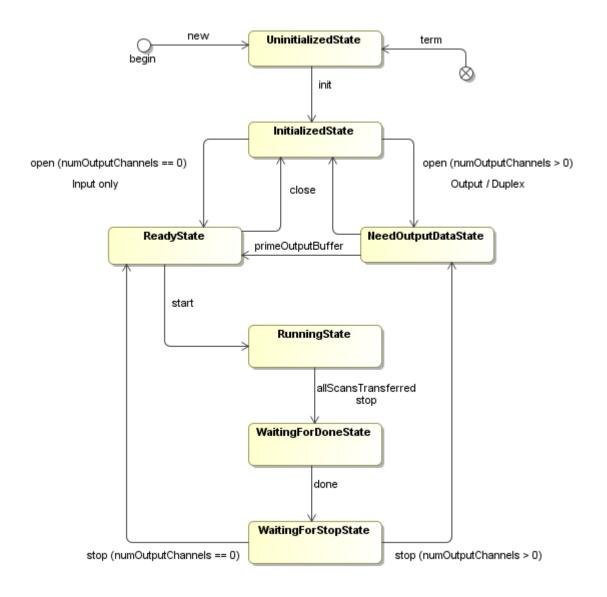

- "State Machine Diagram" on page 5-2
- "Foreground Streaming Sequences" on page 5-5
- "Background Streaming Sequences" on page 5-12

Streaming Sequence Diagrams

These sequence diagrams provide details of timing for streaming functionality. They might be useful for debugging code during development of your adaptor.

•

- "Foreground Streaming Sequences" on page 5-5
- "Background Streaming Sequences" on page 5-12
- "Sequence for Errors and Exceptions" on page 5-21

Foreground Streaming Sequences


This topic includes sequence diagrams for finite analog input and output in the foreground.

In this section...

"Sequence for Finite Foreground Input" on page 5-5 "Sequence for Finite Foreground Output" on page 5-7 "Sequence for Finite Foreground Duplex Channel" on page 5-9


Sequence for Finite Foreground Input

This diagram shows the timing sequence for a finite (fixed-size) analog input in the foreground. It demonstrates the interfaces between a DAQ session, AsyncIO channel, and a DAQ stream when a user is performing finite foreground clocked acquisitions using the session interface.

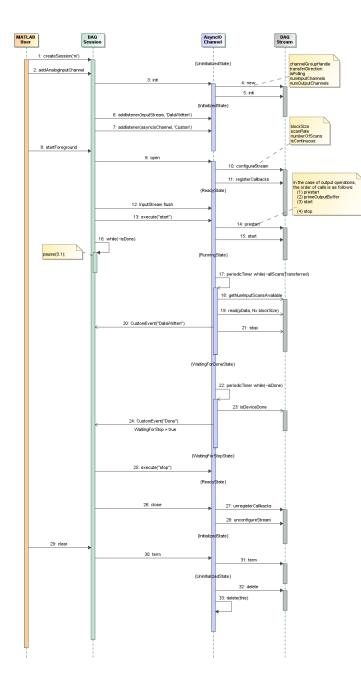


Sequence for Finite Foreground Output

This diagram shows the timing sequence for a finite (fixed-size) analog output in the foreground. It demonstrates the interfaces between a DAQ session, AsyncIO channel, and a DAQ stream when a user is performing finite foreground clocked signal generation using the session interface.

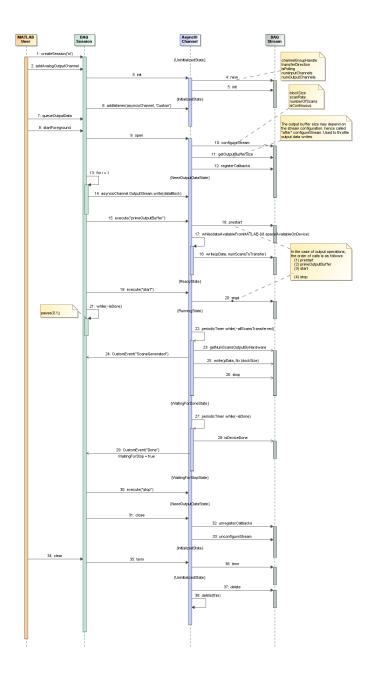

Sequence for Finite Foreground Duplex Channel

This diagram shows the timing sequence for a finite (fixed-size) simultaneous analog input and output in the foreground. It demonstrates the interface between a DAQ session and the AsyncIO channel when a user is performing a finite input/output operation using the session interface

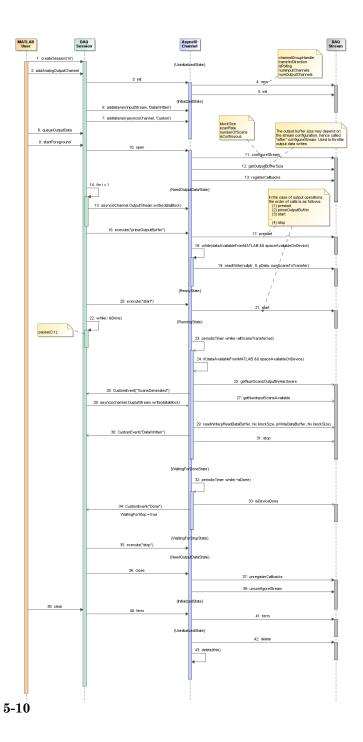

- "State Machine Diagram" on page 5-2
- "Background Streaming Sequences" on page 5-12

Background Streaming Sequences


This topic includes sequence diagrams for analog input and output in the background.

In this section...

"Sequence for Finite Background Input" on page 5-12 "Sequence for Continuous Background Input with Stop" on page 5-14 "Sequence for Finite Background Input with Wait" on page 5-16 "Sequence for Finite Background Input with Stop Race" on page 5-18


Sequence for Finite Background Input

This diagram shows the timing sequence for a finite (fixed-size) analog input in the background. It illustrates the interface between a DAQ session and the AsyncIO channel when a user is performing a finite background clocked acquisition using a the session interface.

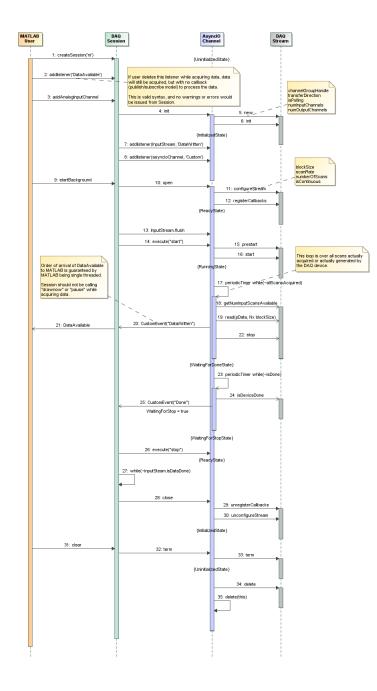


Sequence for Continuous Background Input with Stop

This diagram shows the timing sequence for a continuous analog input in the background, with a stop request while the device is acquiring data. It illustrates the interface between a DAQ session and the AsyncIO channel when a user is performing a continuous background clocked acquisition, then calls stop while data is being acquired.

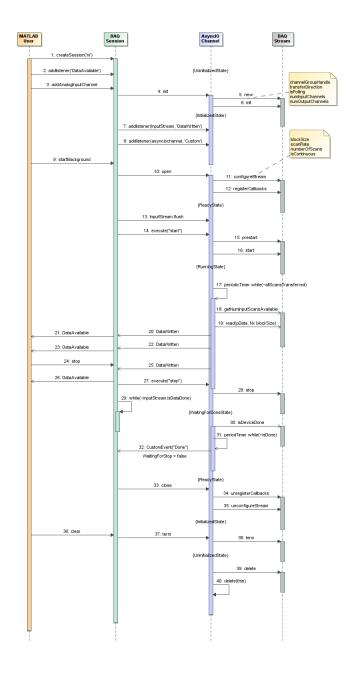


Sequence for Finite Background Input with Wait

This use case revisits the finite background acquisition, when all scans have been acquired and the background operation is naturally stopping, and at the same time, the user issues a stop command while the DAQ AsyncIO plugin is in the "WaitingForDoneState."

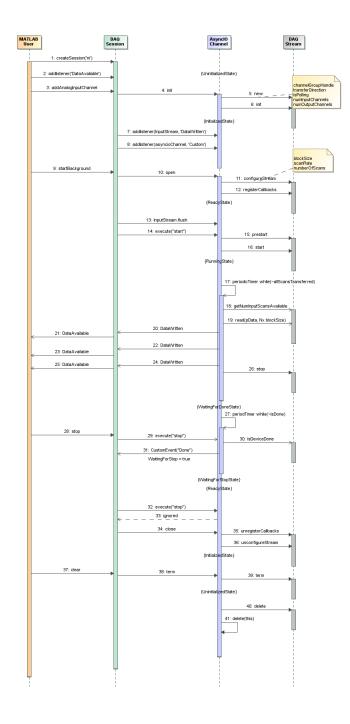

Sequence for Finite Background Input with Stop Race


This situation revisits the finite background acquisition, when all scans have been acquired and the background operation is naturally stopping, and at the same time, the user issues a stop command while the DAQ AsyncIO plugin is in the last iteration of the "RunningState."

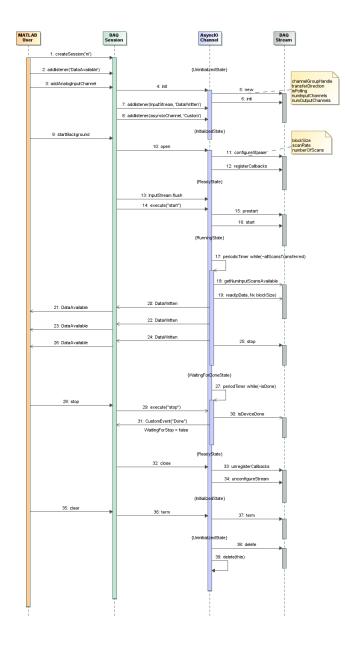

- "State Machine Diagram" on page 5-2
- "Foreground Streaming Sequences" on page 5-5

Sequence for Errors and Exceptions

This sequence diagram summarizes the DataMissed events and error handling.

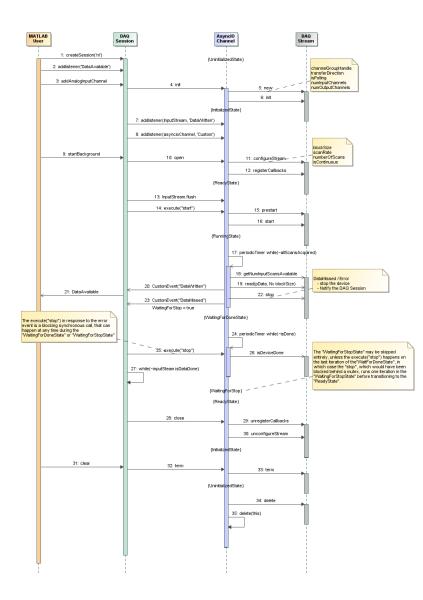

- "Channel Groups" on page 3-13
- "Errors and Exceptions" on page 3-11
- "State Machine Diagram" on page 5-2
- "Foreground Streaming Sequences" on page 5-5
- "Background Streaming Sequences" on page 5-12

Functions — Alphabetical List

buildAdaptor

Build adaptor for third-party data acquisition interface

Syntax

```
daq.sdk.utility.mex.buildAdaptor(adaptorName,customFunc,srcPath,
outputPath)
daq.sdk.utility.mex.buildAdaptor(adaptorName,customFunc,srcPath,
outputPath,vendorLib)
script = daq.sdk.utility.mex.buildAdaptor(____)
```

Description

daq.sdk.utility.mex.buildAdaptor(adaptorName,customFunc,srcPath, outputPath) builds an adaptor for enumerating, configuring, and streaming data to and from a data acquisition device driver.

Note This function requires that your system is configured with Microsoft Visual Studio 2013 or later.

daq.sdk.utility.mex.buildAdaptor(adaptorName,customFunc,srcPath, outputPath,vendorLib) allows you to specify a custom library for the build.

script = daq.sdk.utility.mex.buildAdaptor(____) returns the script used for the build. This can be useful for diagnostic purposes.

Examples

Build Custom Adaptor

Build the custom adaptor named MyAdaptor.

daq.sdk.utility.mex.buildAdaptor('MyAdaptor', 'custom my', 'c:\adaptors\sdk\daqadaptor', 'c:\adaptors\sdk\bin')

Build with Custom Library

Use the custom library MyLibrary for building an adaptor.

```
pathToHeaderAndLib = 'C:\libraries\MyLibrary'
myLibrary.HeaderPath = fullfile(pathToHeaderAndLib,'include');
myLibrary.LibPath = fullfile(pathToHeaderAndLib,'lib');
myLibrary.LibName = 'MyLibrary';
```

buildAdaptor('DemoAdaptor', 'custom_demo', adaptorPath, outputPath, myLibrary);

"Create Your Adaptor from the Demo Adaptor" on page 3-2

Input Arguments

adaptorName - Name of adaptor

char vector | string

Name of the adaptor, specified as a character vector or string.

Example: 'DemoAdaptor'

Data Types: char | string

customFunc — File containing custom functions

char vector or string

Name of the file containing the source code for custom functions, specified as a character vector or string. The file must be in the folder identified by srcPath.

Example: 'custom_demo.cpp'

Data Types: char | string

srcPath — Path to adaptor source folder

char vector or string

Example: 'c:\temp\sdk\daqadaptor'

Data Types: char | string

outputPath — Path to adaptor MEX-file

char vector or string

Path to generated adaptor MEX-file location, specified as a character vector or string.

Example: 'c:\temp\sdk\bin'

Data Types: char | string

vendorLib - Vendor library locations struct

Vendor library locations, specified as a structure containing these three fields:

- HeaderPath a character vector specifying the path to the vendor header.
- LibPat a character vector specifying the path to the vendor static library.
- LibName a character vector specifying the name of the static library, without file extension.

Data Types: struct

Output Arguments

script — Build script
character vector

Build script returned as a character vector. This script indicates what the function ran to build the adaptor.

See Also

Topics

"Create Your Adaptor from the Demo Adaptor" on page 3-2

Introduced in R2017a

enableDemoAdaptorDiscovery

Allow SDK demo adaptor to be enabled for device discovery and usage

Syntax

daq.sdk.utility.enableAdaptorDiscovery

Description

daq.sdk.utility.enableAdaptorDiscovery allows the SDK demo adaptor to be found and used in a data acquisition session. The adaptor is enabled until the end of the MATLAB session or until execution of daqreset.

Examples

Enable the Demo Adaptor

Enable the demo adaptor and view its devices.

```
dagreset;
daq.sdk.utility.enableDemoAdaptorDiscovery;
devices = daq.getDevices
devices =
Data acquisition devices:
index Vendor Device ID
                           Description
1
              MWDevO
                        MathWorks MW314159
      mw
2
      mw
             MWDev1
                        MathWorks MW314159
                        MathWorks MW628318
3
      mw
             MWDev2
```

- "Demo Adaptor Description" on page 2-2
- "Enable the Demo Adaptor" on page 2-5
- "Session Workflows with the Demo Adaptor" on page 2-6

• "Create Your Adaptor from the Demo Adaptor" on page 3-2

See Also

See Also

Functions

daqreset

Topics

"Demo Adaptor Description" on page 2-2 "Enable the Demo Adaptor" on page 2-5 "Session Workflows with the Demo Adaptor" on page 2-6 "Create Your Adaptor from the Demo Adaptor" on page 3-2

Introduced in R2017a